
## ALCOSENS SYSTEMS ONLINE DETERMINATION OF ALCOHOL

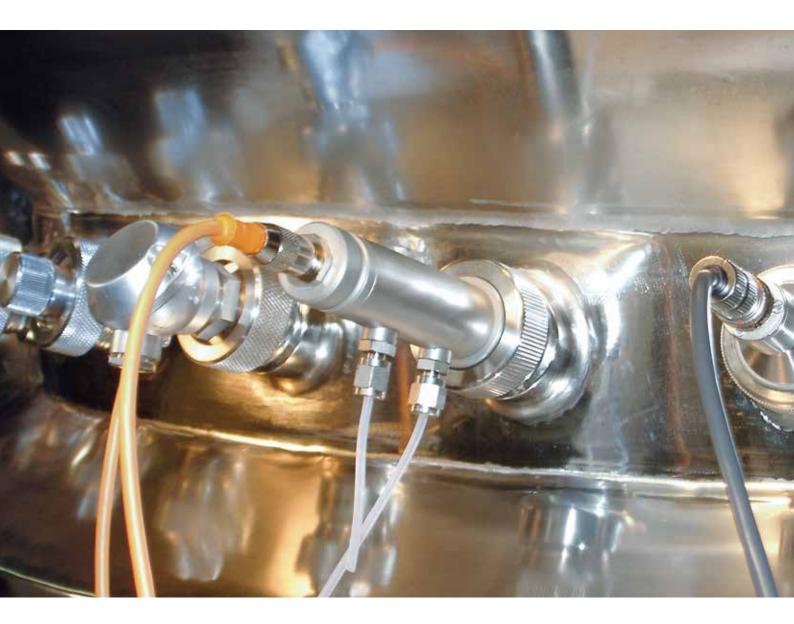

ALCOSENS PROBE AND TRANSMITTER



## **Characteristics**:

- Online determination of alcohol or other volatiles
- Measuring of alcohol concentration in liquids
- Design for cleaning (CIP) and sterilisation (SIP)
- Customized probe geometries available
- Transmitter/controller for up to 8 probes
- Touchscreen for visualisation and calibration
- Interface to central PLC and SCADA systems
- GMP documentation on demand

# ALCOSENS SYSTEMS ETHANOL AND METHANOL MONITORING AND CONTROL

### **ALCOSENS** Applications

ALCOSENS systems are the standard automation equipments for alcohol determination in a number of fermentations such as industrial yeast fermentation and acetic acid fermentation. The ALCOSENS technology is also widely used for the online measurement of methanol and the sophisticated control of processes with recombinant methylotrophic yeasts.

- Methanol monitoring and control in *Pichia pastoris* fermentation in fed batch mode
- Ethanol control in baker's yeast production
- Control of ethanol concentration in acetic acid fermentation
- Process optimization in brewers yeast propagation
- Observing dealcoholization of beer
- Applications in bioethanol plants



Transmitter S7 Control Cabinet with Touch Panel

### **Probe Geometries and Transmitter Configurations**

ALCOSENS probes are available in different geometrical configurations and with different port adapters. In combination with the TRANSMITTER S7 MULTI for up to 8 probes the system can be used for small scale parallel bioreactors, too.

#### **General Specifications**

Measuring range:	0.06-6% (v/v)
Transmitter output:	4-20 mA / 0 - 10 V
Sensor:	Metal oxide semiconductor
Probe body:	AISI 316L
Membrane:	Silicone
Probe adaptor:	19 mm / 25 mm port
Lengths of probe:	105/120/220/320mm
Max. pressure:	2.5 bar
Temperature:	20 – 143 °C (for sterilisation only)
Carrier gas supply:	20 mL / min (synthetic or purified air)



ALCOSENS Standard Probe for 25 mm Port, side wall mounting



ALCOSENS Customized Probe for 19 mm Port, headplate mounting



ALCOSENS G Customized Probe for GMP Applications

Frings	Ferm Ferm Ferm	enter 2 6.00 enter 3 6.00 enter 4 2.15 enter 5 3.34	
Calibration		Settings	System
Trend	re	Trend	Alarm Message
9		nperature ( *C )	

03,05,2012 10:29:3	7 Fermenter #6
Oter: Vef %	1 12 12 5 1 100 2200 5 1 100 5 100
(t) Qh	30.0 < Onive 1 (100 1 12)
Pecalibration :	a chen a co mi fa
Ohm NC : Ohm NC : Online Yol % :	Ofm Ofm Ofm Vol %
di sorten	ACI

Interface of ALCOSENS TRANSMITTER S7 MULTI with Overview, Calibration and Trend Window

Heinrich Frings GmbH & Co. KG Boschstraße 32, D-53359 Rheinbach Phone: +49 (0)22 26 89 29 - 400 Fax: +49 (0)22 26 89 29 - 401 marketing@frings.com www.frings.com